


# ColorFlex® EZ Coffee Quick Start Guide

A60-1016-265 ver 1.1

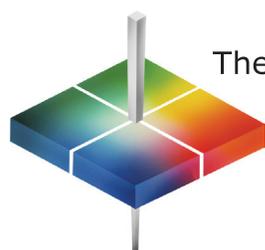

The world's true measure of color

# HunterLab

## A SETTING UP THE INSTRUMENT

1. Connect the AC adapter to the left port on the back of the ColorFlex EZ.



2. Plug the power cord into a wall outlet.



3. If a printer or other USB device is to be used, plug into one of the two USB ports on the side of the ColorFlex EZ.



4. Turn on the ColorFlex EZ by pressing **GO** (⚡) on the button pad.



Now your instrument is ready for Standardization.

5. (Optional) Using HunterLab's EasyMatch QC software to control ColorFlex EZ

Connect the USB cable from the rear Mini-B plugin on the ColorFlex EZ to a standard USB plugin on the PC. Load EasyMatch QC on the computer and connect to the ColorFlex EZ as a sensor.

## B STANDARDIZATION

1. Press the **DOWN** arrow to select **STANDARDIZE** from the **MAIN MENU**. Press **GO** to select.



2. Open the standards box and place the black glass at the port. Press **GO**.



3. Place the calibrated instrument white tile at the port. Press **GO**. You are now ready to read samples.



Now your instrument is ready to Read Samples.

## C READING SAMPLES

1. Select **READ** from the **MAIN MENU** and then press **GO**.



2. Press the **UP/DOWN** arrow to select a setup from the list. Press **GO** to continue.



3. Prepare your sample and center it at the port.



4. Press **GO** to read and the color measurement will be displayed.



To view the next Coffee score, press the UP arrow.

5. Press **DOWN** arrow to **SAVE** the reading to memory and **PRINT** the reading to a USB printer.



For more detailed instruction on using the instrument, please refer to the ColorFlex EZ Coffee User's Manual.

Visit [www.hunterlab.com](http://www.hunterlab.com) for information on all our products.